
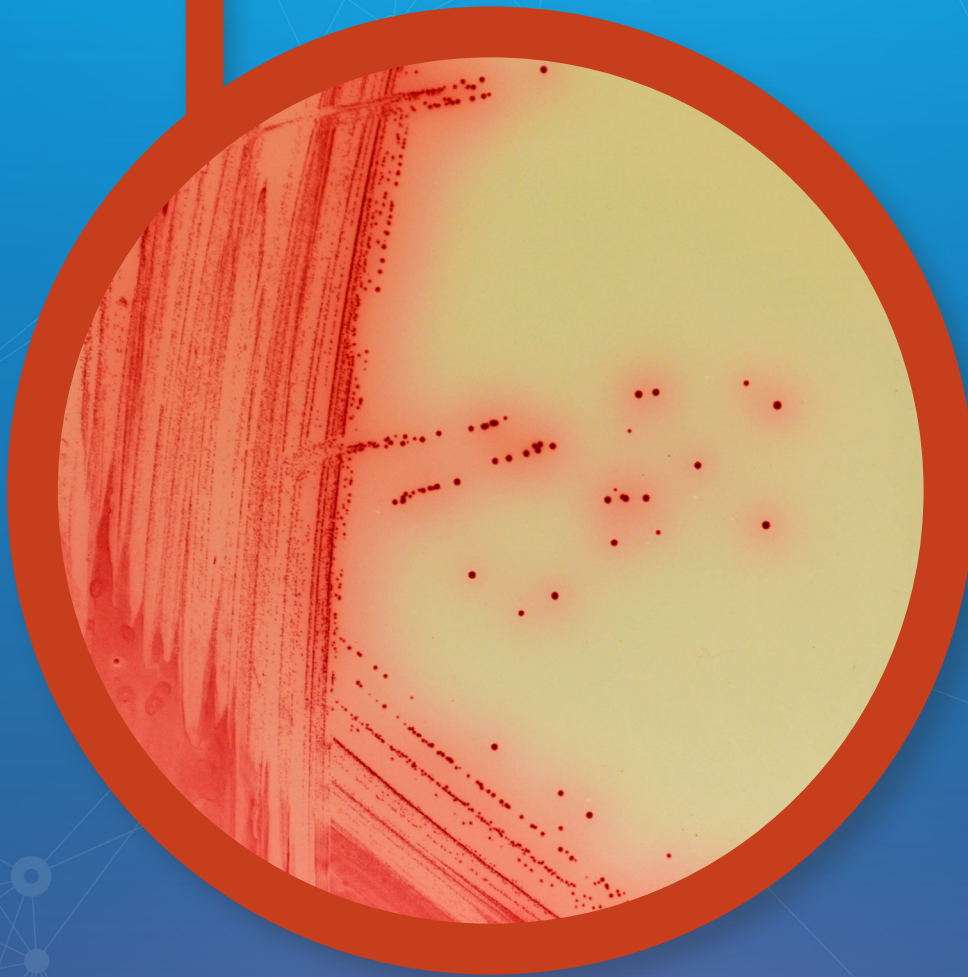
# CHROMagar™ Mycoplasma

Click below:
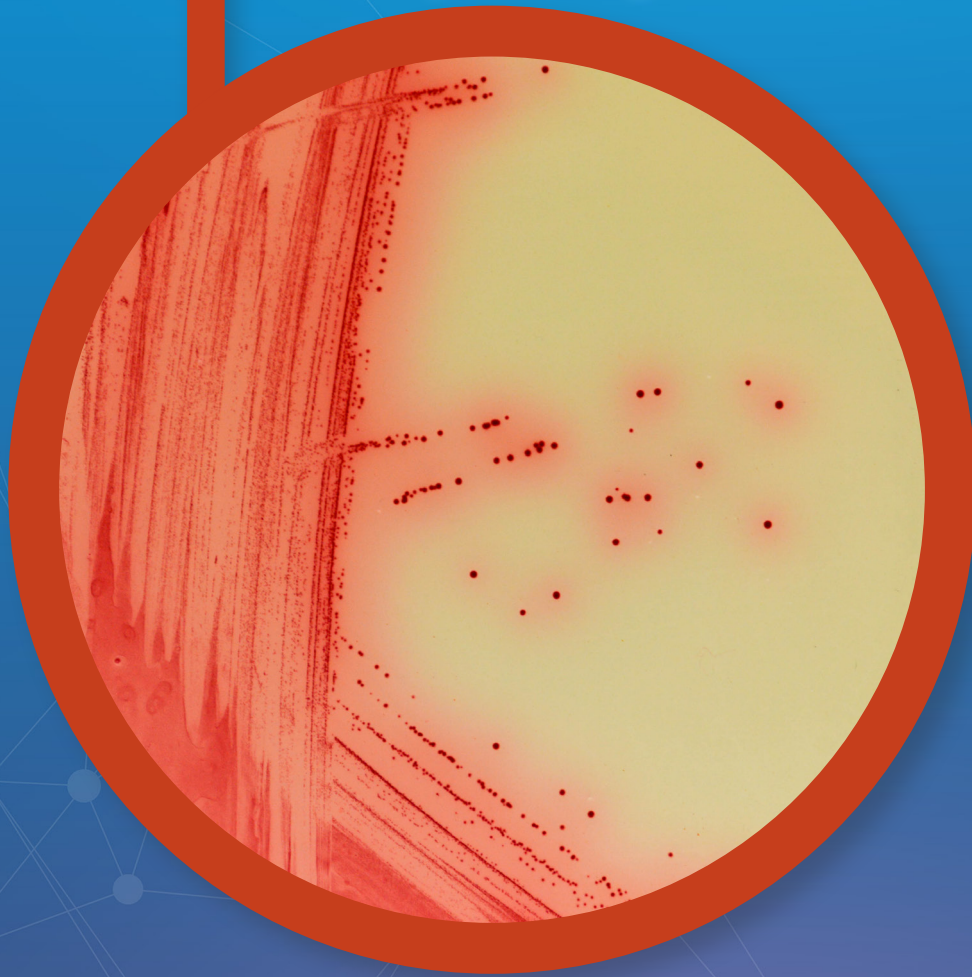
# ● CHROMagar™ Mycoplasma

For detection of *Mycoplasma bovis*

## For detection of *Mycoplasma bovis*

### Intended Use

Bovine infections can result in a variety of clinical signs, such as pneumonia mainly but also arthritis, mastitis and keratoconjunctivitis, none of which are *M. bovis*-specific. Laboratory diagnosis is therefore important.

Identification of *M. bovis* from nose swabs, broncho-alveolar lavage samples can be done by streaking the sample directly onto CHROMagar™ Mycoplasma agar plates.

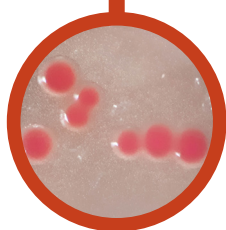
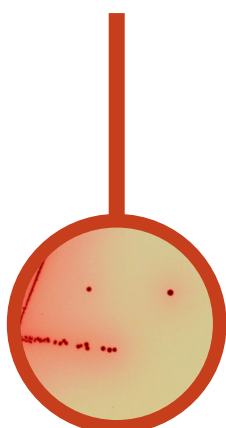
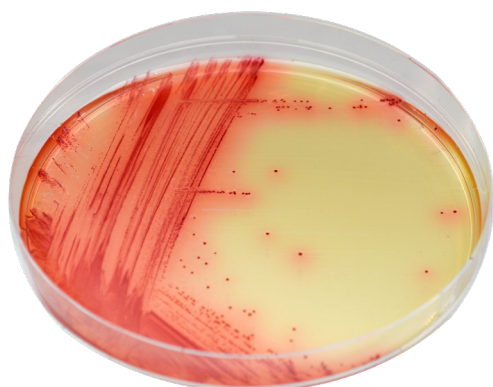
The detection of *M. bovis* by red coloration allows an easy detection with a fried egg aspect under binocular loupe after 3-7 days of incubation at 37 °C under CO<sub>2</sub> atmosphere.

### Medium Performance

- 1 **EASY TO READ AND INTERPRET**  
With CHROMagar™ Mycoplasma, *Mycoplasma* colonies have an intense red coloration and a fried egg appearance - making detection very easy.
- 2 **DETECTION**  
Detection is possible between 3 to 7 days under CO<sub>2</sub> atmosphere.
- 3 **SIMPLE METHOD**  
Identification of *M. bovis* from nasal swabs and broncho-alveolar lavage samples can be performed by plating the sample directly onto the plates.

### Medium Description

<b>Powder Base</b>	Total .....	43.8 g/L
	Agar .....	10.0
<b>Supplement (S1)</b> included in this reference	Peptone and yeast extract .....	25.0
	Mineral Salts .....	6.3
<b>Supplement (S2)</b> included in this reference	Growth factors .....	2.2
	Chromogenic and selective mix .....	0.3
<b>Supplement (S3)</b> included in this reference	Storage at 15/30 °C - pH: 7.6 +/- 0.2	
	Shelf Life .....	> 24 months
Usual Samples	Powder .....	1 g/L
	Storage at 15/30 °C	
Procedure	Shelf Life .....	> 24 months
	Powder .....	1 g/L
Scientific Publications on this product: available on <a href="http://www.CHROMagar.com">www.CHROMagar.com</a>	Storage at 2/8 °C	
	Shelf Life .....	> 18 months
Please read carefully the instructions for use (IFU document) available on : <a href="http://www.CHROMagar.com">www.CHROMagar.com</a>		

### Plate Reading

- *Mycoplasma bovis*  
→ red
- *Mycoplasma bovis*  
under binocular loupe  
→ red with a fried egg  
morphology

### Order References

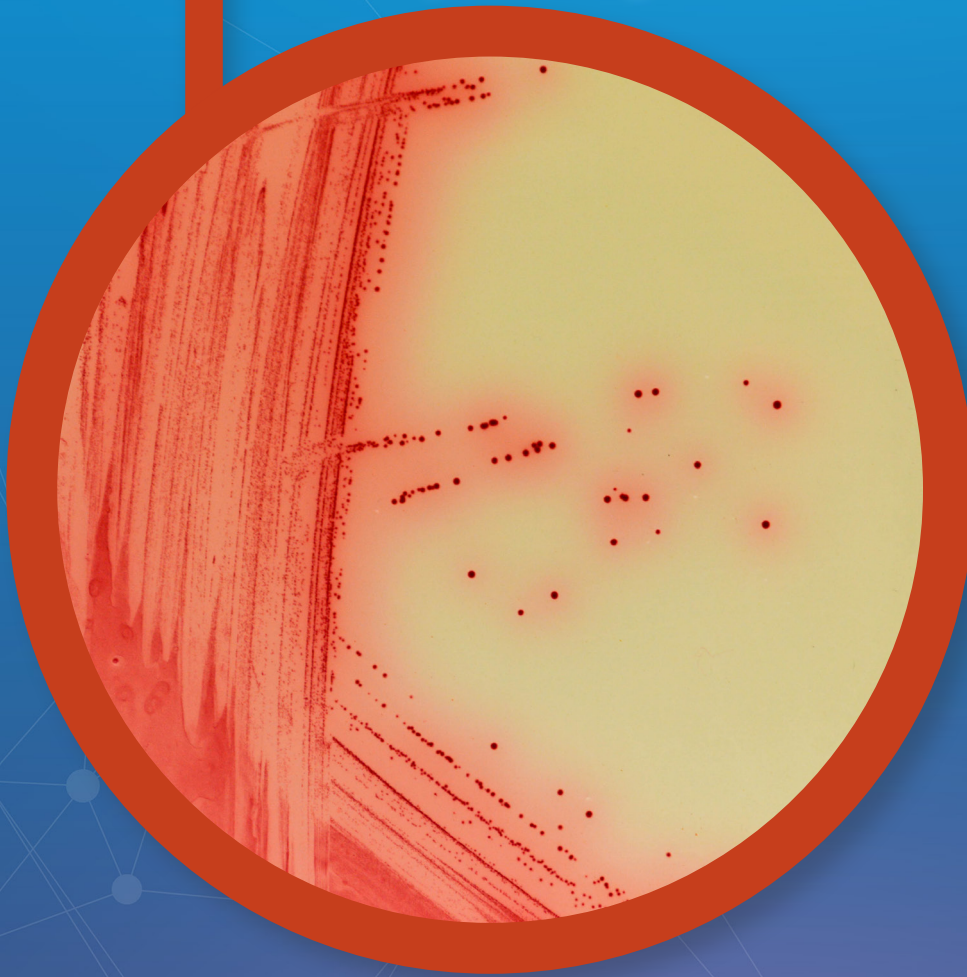
Please use these product references when contacting your local distributor:

1000 mL pack ..... MB640

(Included in this reference: powder base MB640(B) + supplement MB640(S1) + supplement MB640(S2) + supplement MB640(S3))

Manufacturer: CHROMagar  
4 place du 18 juin 1940 75006 Paris - France  
Email: [CHROMagar@CHROMagar.com](mailto:CHROMagar@CHROMagar.com)  
Website: [www.CHROMagar.com](http://www.CHROMagar.com)  
Find your nearest distributor on  
[www.CHROMagar.com/contact](http://www.CHROMagar.com/contact)

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